

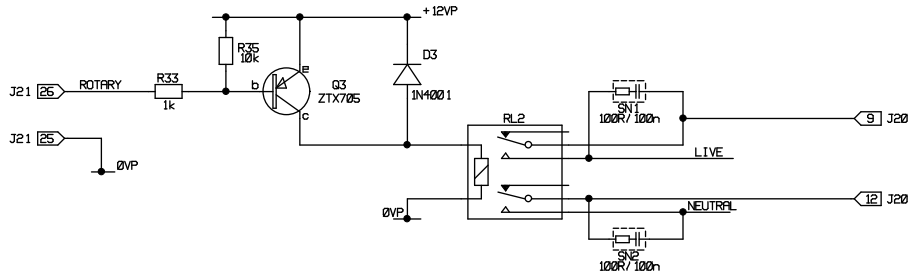
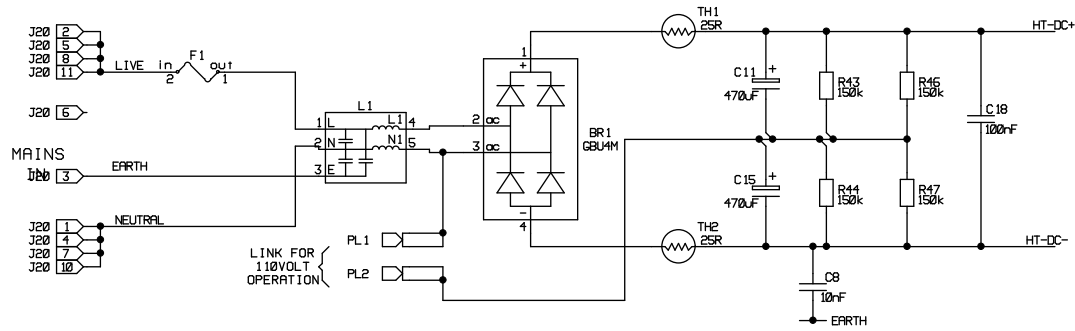


MODIFICATIONS		EMITECH	
TITLE: Standard H.T. Supply PCB		ISSUE: 2	
DRAWN DJR	DATE: 07/04/2000	DRG NO. A3/001/210	SHEET 1 OF 2

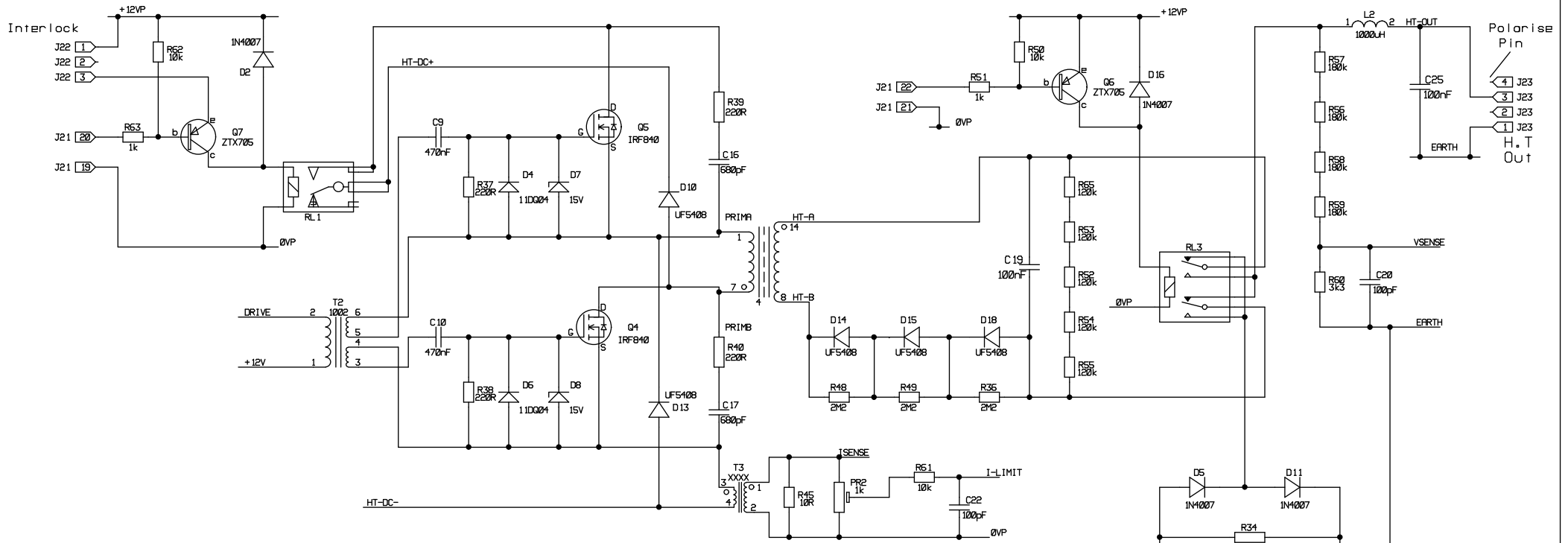



MODIFICATIONS		EMITECH	
TITLE: Standard H.T. Supply PCB		ISSUE: 2	
DRAWN DJR	DATE: 07/04/2000	DRG NO. A3/001/210	SHEET 2 OF 2

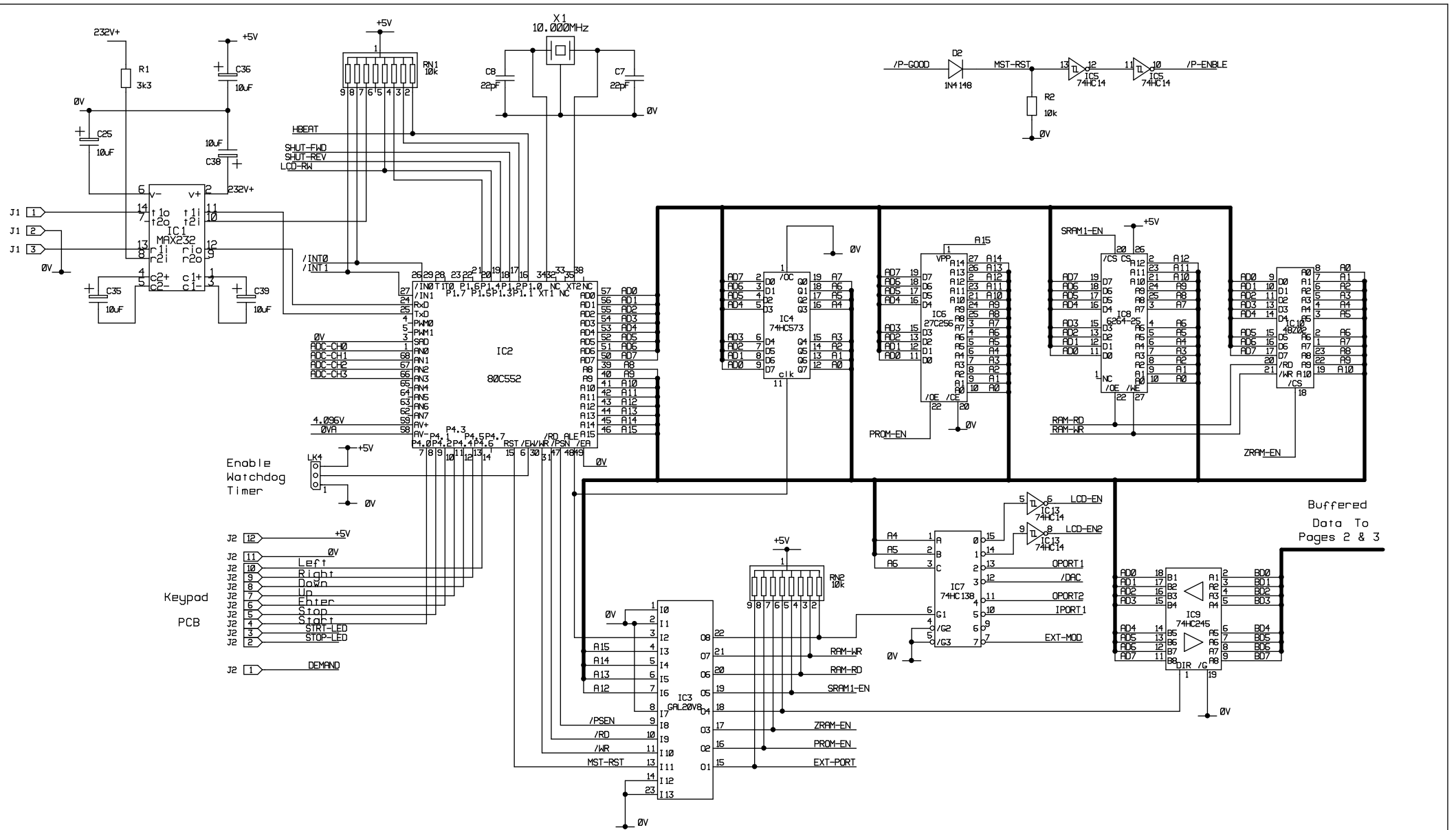








Buffered
Data To
Pages 2 & 3

MODIFICATIONS		EMITECH	
TITLE:		K550X Control PCB Circuit Schematic	
DRAWN DJR		DATE: 15/02/2000	DRG NO. A3/107/001
		ISSUE: 1	SHEET 1 OF 4







Buffered I/O Data Bus

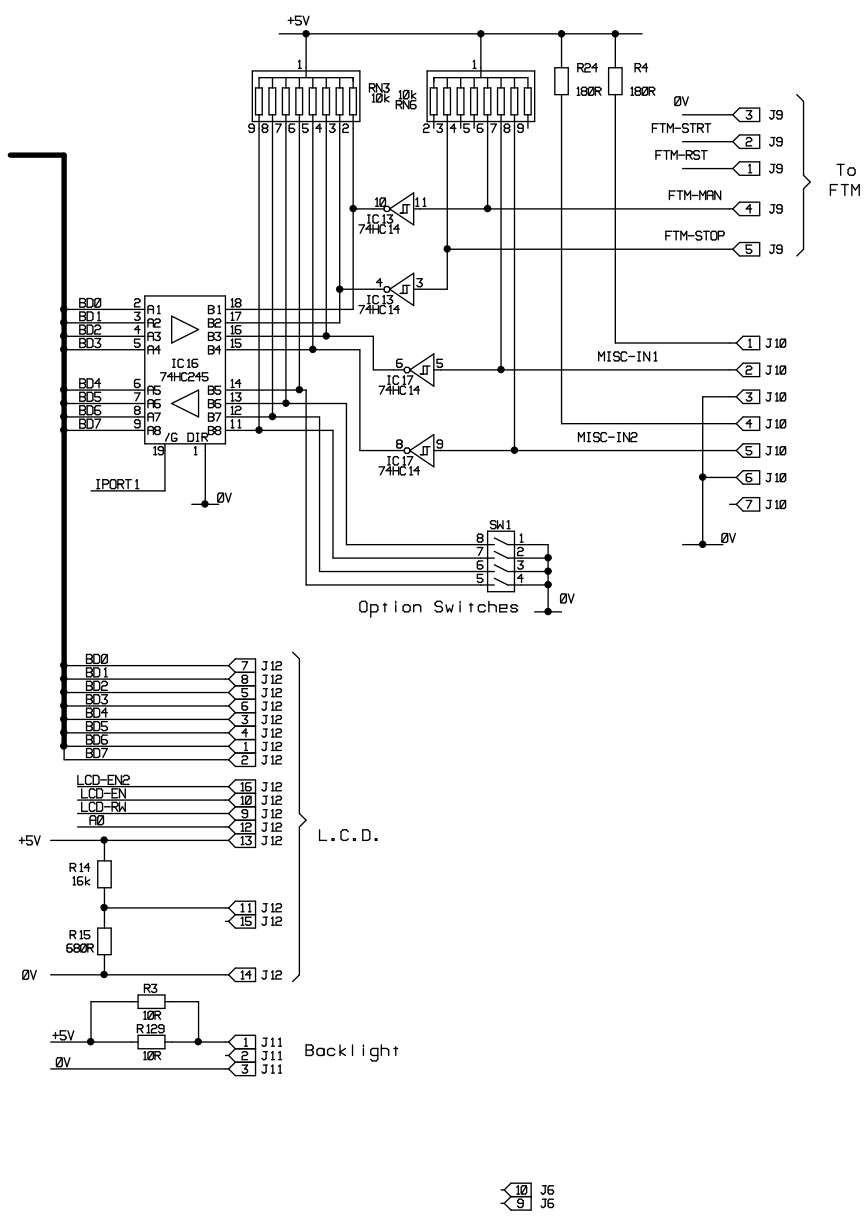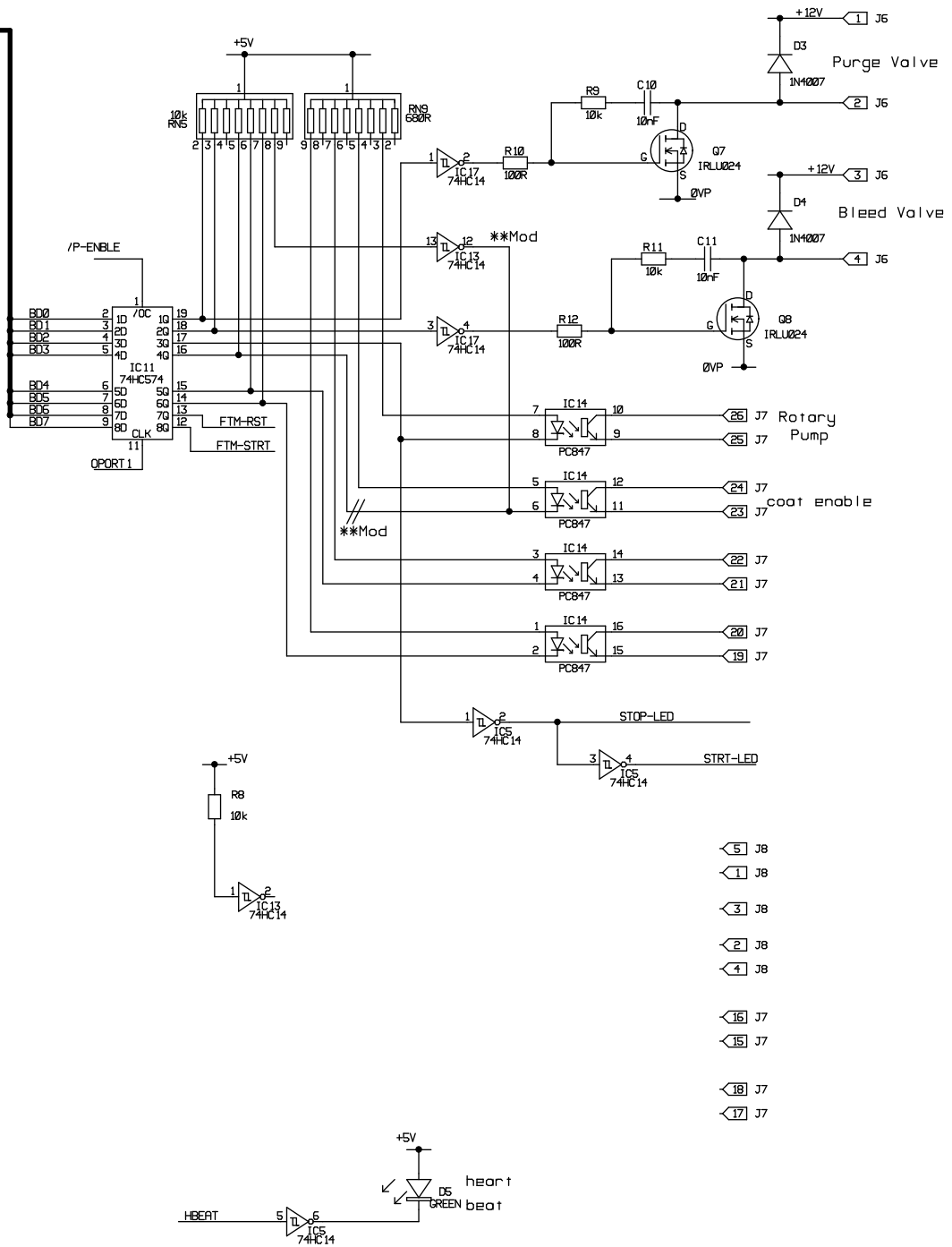


MODIFICATIONS		EMITECH	
TITLE:		K550X Control PCB Circuit Schematic	
DRAWN DJR		DATE: 15/02/2000	DRG NO. A3/107/001
		ISSUE: 1	SHEET 2 OF 4

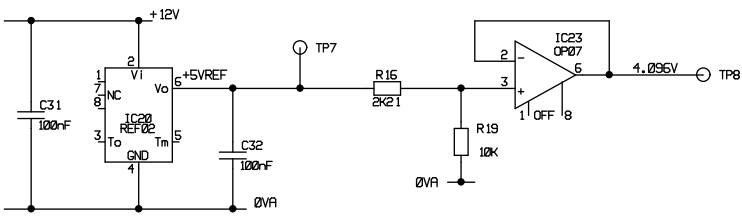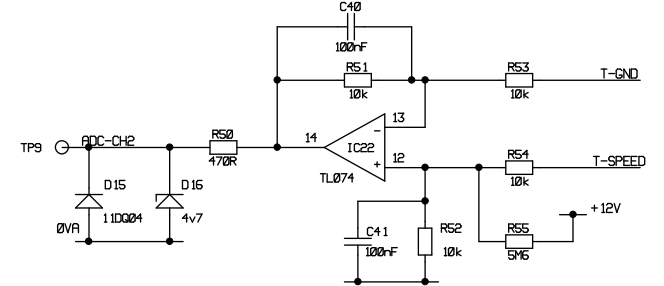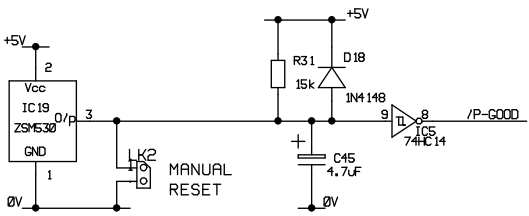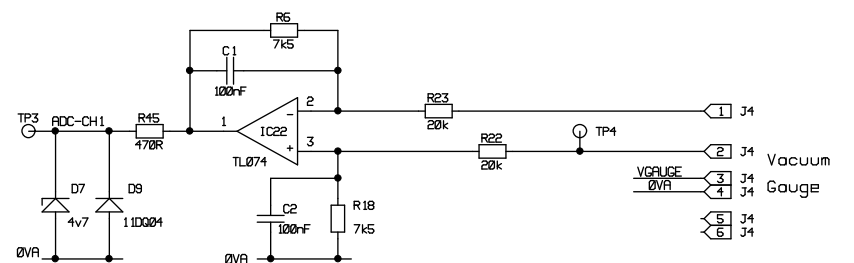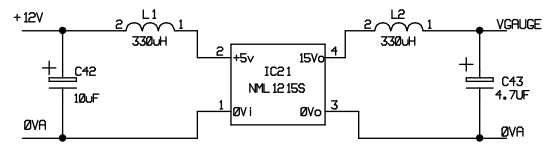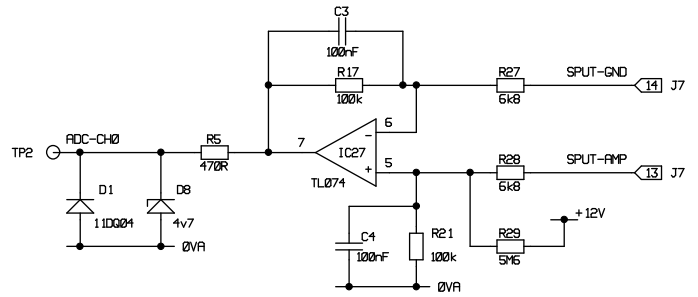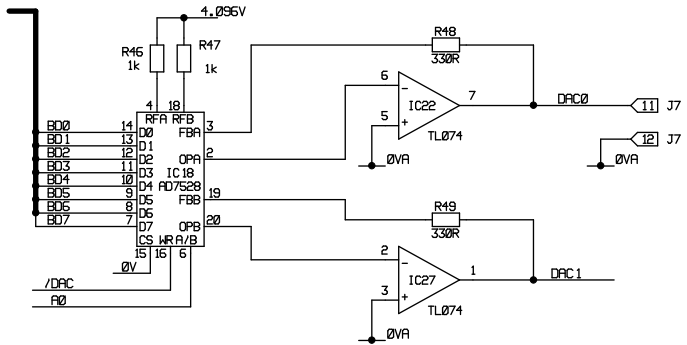

MODIFICATIONS		EMITECH	
TITLE:		K550X Control PCB Circuit Schematic	
DRAWN DJR		DATE: 15/02/2000	DRG NO. A3/107/001
		SHEET 3 OF 4	

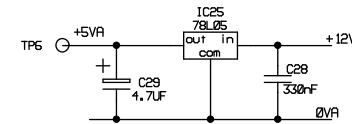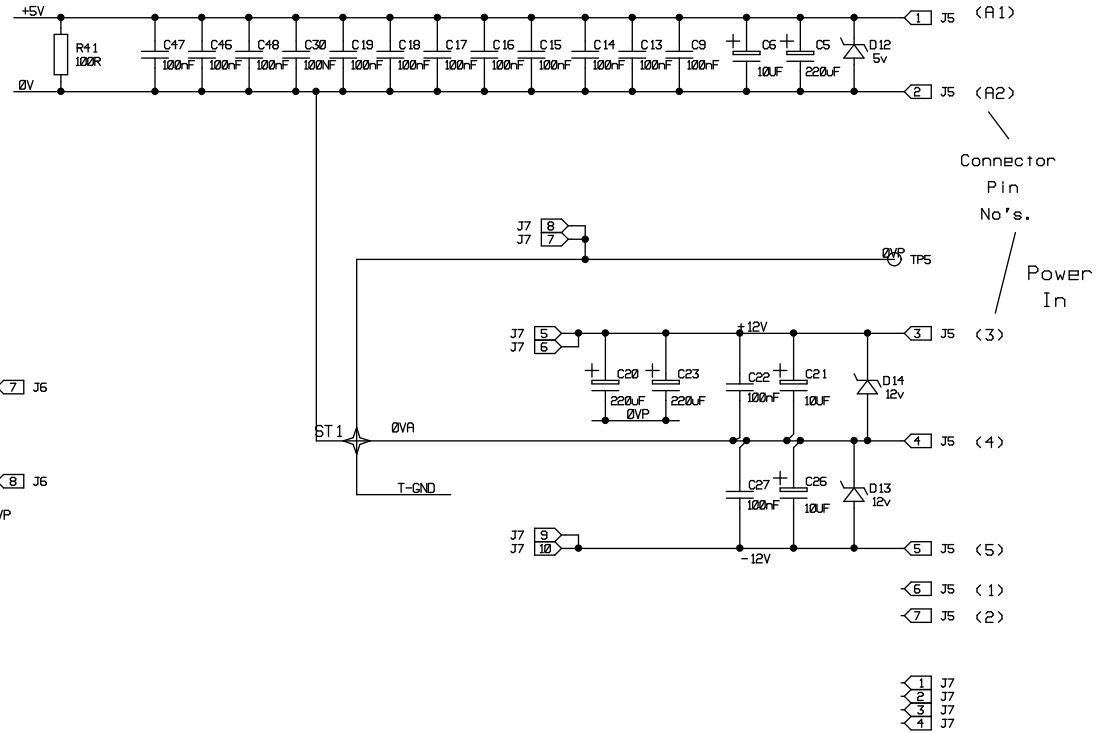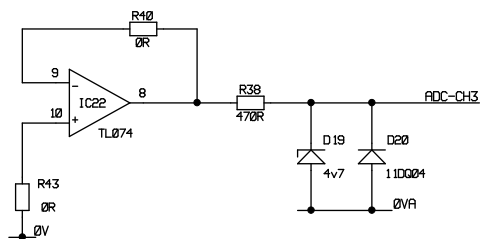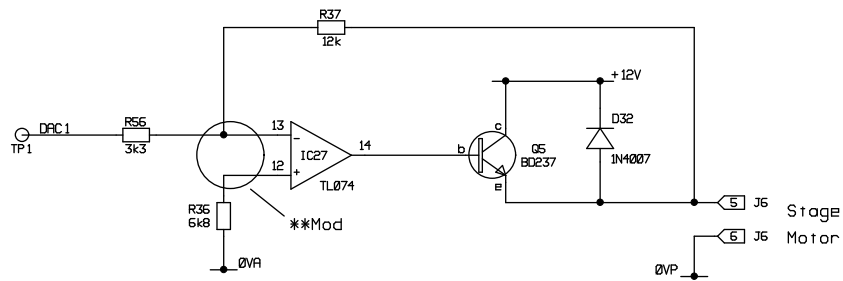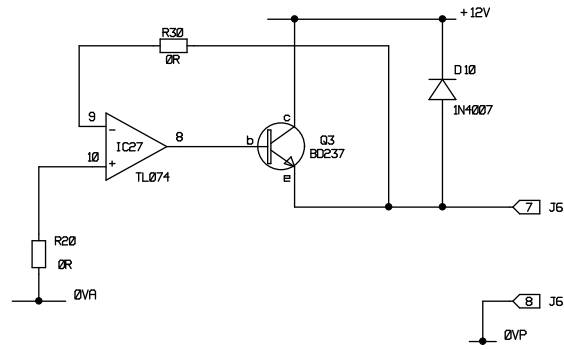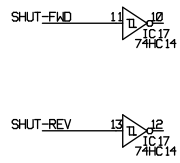

MODIFICATIONS		EMITECH	
TITLE:		K550X Control PCB Circuit Schematic	ISSUE: 1
DRAWN DJR	DATE: 15/02/2000	DRG NO. A3/107/001	SHEET 4 OF 4

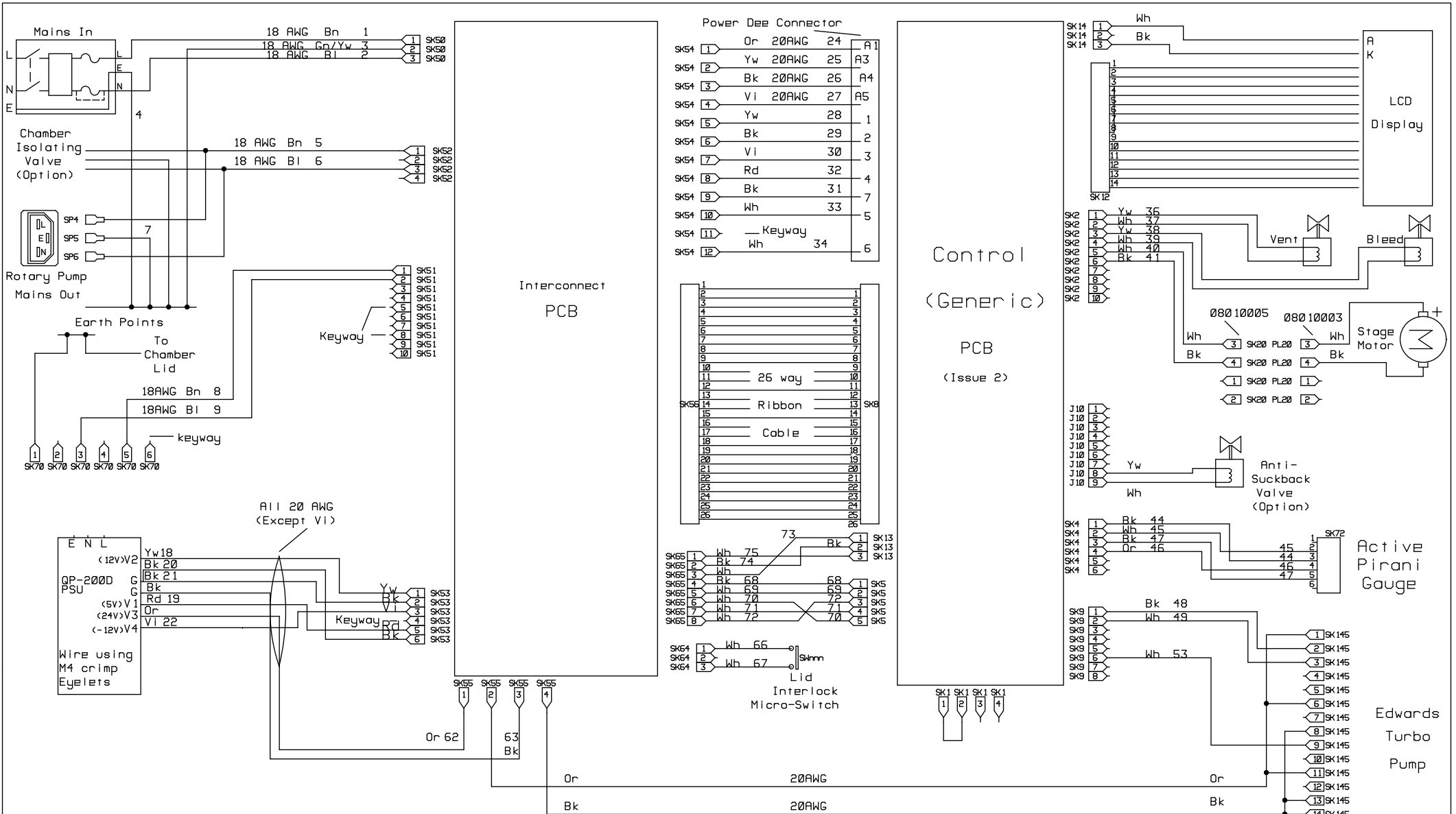


All Wires 24AWG unless otherwise specified


Colour code	
Yw	- Yellow
Wh	- White
Bl	- Blue
Bk	- Black
Br	- Brown
GrY	- Green/Yellow
Gr	- Grey

MODIFICATIONS Ref CAR 0527, 0532		EMITECH	
TITLE: K650XT Wiring Schematic		ISSUE: 2	
DRAWN CJH	DATE: 19/5/2006	DRG NO. A3/307/016	SHEET 1 OF 2

Sputter
Supply
PCB

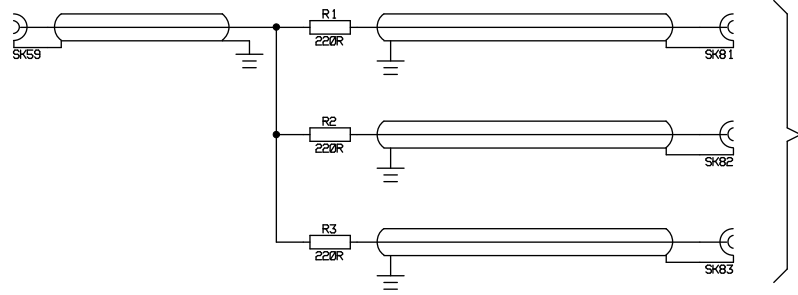

Sputter
Target
Heads

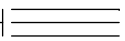MODIFICATIONS		EMITECH 	
TITLE: K650XT Wiring Schematic		ISSUE: 2	
DRAWN CJH	DATE: 19/5/2006	DRG NO. A3/307/016	SHEET 2 OF 2