



Quorum Technologies

SC7610 Sputter Coater

Operating Manual



Document Number OM-SC7610

Issue 1 (01/02)





Quorum Technologies

Quorum Technologies Ltd is the owner and manufacture of the preparation equipment.



range of EM

Unit 15A
Euro Business Park
New Road
Newhaven
East Sussex
U.K.
BN9 0DQ
Email :

tel: ++44(0) 1273 510535
fax: ++44(0) 1273 510536

sales@quorumtech.com
<http://www.quorumtech.com>

For further information regarding any of the other products designed and manufactured by Quorum Technologies, contact your local representative or directly to Quorum Technologies at the address above.

- Carbon and Sputter Coaters
- Plasma Reactor for ashing and etching
- High Vacuum Bench Top Evaporators
- Cryo Transfer Systems
- Critical Point Dryers
- Service and Spares

Disclaimer

The components and packages described in this document are mutually compatible and guaranteed to meet or exceed the published performance specifications. No performance guarantees, however, can be given in circumstances where these component packages are used in conjunction with equipment supplied by companies other than Quorum Technologies



1 Contents

1.1 Manual Layout

This Operating Manual is divided up into the following major section, each chapter dealing with specific topics, as follows:

Chapter 1 - Contents

Chapter 2 - Health and Safety

General section which applies to all Quorum Technologies products detailing the very important issues of Health and Safety applicable when using sample preparation equipment.

Chapter 3 - Introduction

Introduces this manual.

Chapter 4 - General Description

Identifies each of the equipment items and provides an overview of their functions and how they work.

Chapter 5 - Installation

Instructions on how this Instrument should be installed and the connections which should be made between the equipment items.

Chapter 6 - Operation

Instructions on how to start-up and run the instrument.

Chapter 7 - Maintenance

Instructions on how to check the system is functioning correctly, and how to change consumable items. Details of appropriate Spare parts.

Chapter 8 - Fault Finding

Information on how to identify faults in the system and how to rectify these faults.

Chapter 9 - Agents

List of agents supporting Quorum Technologies products.

Chapter 10 - Index

1.2 Chapter Contents

Chapter 1 - Contents	3
1.1 Manual Layout	3
1.2 Chapter Contents	4
1.3 Illustrations	6
1.4 Tables	6
Chapter 2 - Health and Safety	7
2.1 Control of Substances Hazardous to Health (COSHH)	7
2.2 Safety Policy	7
2.3 Conformity	7
2.4 Servicing	8
2.4.1 Disclaimer	8
2.4.2 Operators and Service Engineers	8
2.5 Hazard Signals and Signs	8
2.5.1 Hazard Signal Words	8
2.5.2 Hazard Labels used on Equipment	8
2.5.3 Hazard Warning Labels used in Equipment Manuals	9
2.5.4 Instrument Functionality Signs	9
2.5.5 Serious Damage to Instruments	9
2.5.6 Hazard to Operator	9
2.6 Risk Analysis	10
2.6.1 Personal Operational Risks	10
2.6.2 Hazardous Materials	10
2.7 Good Working Practices	11
2.8 SC7610 Sputter Coater Specific Safety Hazards	11
2.8.1 Contamination	11
Chapter 3 - Introduction.....	12
3.1 Return of Goods.....	12
3.2 Returns Procedure.....	12
Chapter 4 - Description	13
4.1 Equipment	13
4.1.1 Optional Items	13
4.1.2 Accessories	13
4.2 Overview	14
4.3 Technical Specification	14
4.3.1 SC7610 Sputter Coater Specification	14
4.4 Physical Description	15
4.4.1 Operational Description	16
4.4.2 Pumping Requirements	16
4.5 Interlocks	16
4.6 SC7610 Panel Details	17
4.6.1 SC7610, Front Panel Controls	17
4.6.2 SC7610, Rear Panel Connections	18
Chapter 5 - Installation	19
5.1 Unpacking Checklist	19
5.1.1 Preparation	19
5.2 SC7610 Installation	20
5.2.1 Connections	20

Chapter 6 - Operation	22
6.1 Test Procedure	22
6.2 Coating Process	23
6.3 Coating Specimens	23
6.4 Sputtering Rates	25
Chapter 7 - Maintenance	26
7.1 Maintenance - General	26
7.2 Cleaning	26
7.2.1 Vacuum Chamber Cleaning	26
7.3 Target Replacement	27
7.3.1 Target Removal	27
7.3.2 Refit New Target	27
7.4 No Discharge Current	28
7.5 Spare Parts	28
Chapter 8 - Fault Finding	29
8.1 Trouble Shooting / Fault Finding	29
8.2 Fault Prevention	29
Chapter 9 - Agents	30
Chapter 10 - Index	31

1.3 Illustrations

Figure 2.1 - Hazard Warning Symbols 8
Figure 2.2 - International Warning Symbols 9
Figure 2.3 - Typical Signs as shown in this Manual 9
Figure 2.4 - Typical Signs as shown in this Manual 9
Figure 2.5 - Typical Warnings, as shown in this Manual 9
Figure 4.1 - SC7610 Sputter Coater 15
Figure 4.2 - SC7610 Sputter Coater Operation 16
Figure 4.3 - SC7610 Sputter Coater, Front Panel Controls 17
Figure 4.4 - SC7610 Sputter Coater, Rear Panel Connections 18
Figure 5.1 - SC7610 Connections 21
Figure 6.1 - SC7610 Sputtering Rates 25
Figure 7.1 - Target Replacement 27

1.4 Tables

Table 2.1 - Personal Operational Risks 10
Table 4.1 - SC7610 Sputter Coater, Front Panel Control Descriptions 17
Table 4.2 - SC7610 Sputter Coater, Rear Panel Connection Descriptions 18
Table 7.1 - Spare Parts 28
Table 8.1 - Trouble Shooting 29
Table 8.2 - Fault Prevention 29

2 Health and Safety

Safety is very important when using any instrumentation and this chapter should be read by all users of our equipment.

This section of the Manual applies to all surface analysis and surface preparation equipment supplied by Quorum Technologies Polaron range of products, not just the particular instrument for which the manual refers.

Included in this chapter are details on warning notations, good working practices and information on European Community (EC) legislation regarding **“Control Of Substances Hazardous to Health”** (COSHH) and risk analysis.

2.1 Control of Substances Hazardous to Health (COSHH)

The E.C. legislation regarding the “Control of Substances Hazardous to Health” requires Quorum Technologies to monitor and assess every substance entering or leaving their premises. Consequently any returned goods of whatever nature must be accompanied by a declaration form available from Quorum Technologies, reference number SP-100. Without this declaration Quorum Technologies reserves the right not to handle the substance/item. Also in accordance with E.C. regulations we will supply on request hazard data sheets for substances used in our instruments.

2.2 Safety Policy

This section contains important information relating to all health and safety aspects of the equipment. As such it should be read, and understood, by all personnel using the instrument whether as an operator or in a service capacity.

Quorum Technologies is committed to providing a safe working environment for its employees and those that use it's equipment and conducts its business responsibly, and in a manner designed to protect the health and safety of its customers, employees and the public at large. It also seeks to minimise any adverse effects that its activities may have on the environment.

Quorum Technologies regularly reviews its operations to make environmental, health and safety improvements in line with UK and European Community legislation.

The equipment has been designed as free-standing bench mounted instruments. Quorum Technologies cannot be held responsible for any damage, injury or consequential loss arising from the use of its equipment for any other purposes, or any unauthorised modifications made to the equipment.

All service work carried out on the equipment should only be undertaken by suitably qualified personnel. Quorum Technologies is not liable for any damage, injury or consequential loss resulting from servicing by unqualified personnel. Quorum Technologies will also not be liable for damage, injury or consequential loss resulting from incorrect operation of the instrument or modification of the instrument.

2.3 Conformity

This instrument is supplied in a form that complies with the protection requirements of the EC Electromagnetic Compatibility Directive **89/336/EEC** and the essential health and safety requirements of the low voltage directive **72/23/EEC** both as amended by **92/31/EEC**. Any modifications to the equipment, including electronics or cable layout may affect the compliance with these directives.

2.4 Servicing

2.4.1 Disclaimer

All service work on the equipment should be carried out by qualified personnel. Quorum Technologies cannot be liable for damage, injury or consequential loss resulting from servicing from unqualified personnel. Quorum Technologies will also not be liable for damage, injury or consequential loss resulting from incorrect operation of the instrument or modification of the instrument.

2.4.2 Operators and Service Engineers

A normal operator of the equipment will not be trained in or qualified for service work on the equipment and may cause a hazard to himself/herself or others if such work is attempted. Operators should therefore restrict themselves to the normal operation of the equipment and not by removing covers from the electronic equipment or dismantling of the instruments.

Service Engineers who are suitably trained to assess and isolate electrical, mechanical and vacuum hazards should be the only personnel who access the equipment.

2.5 Hazard Signals and Signs

2.5.1 Hazard Signal Words

The standard three hazard signal words are defined as follows:

- ◆ **DANGER** - *imminently* hazardous situation or unsafe practice that, if not avoided, *will* result in death or severe injury.
- ◆ **WARNING** - *potentially* hazardous situation or unsafe practice that, if not avoided, *could* result in death or severe injury.
- ◆ **CAUTION** - *potentially* hazardous situation or unsafe practice that, if not avoided, *may* result in minor or moderate injury or damage to equipment.

2.5.2 Hazard Labels used on Equipment

Several hazard symbols may be found on the equipment, they are shown below with their meaning:

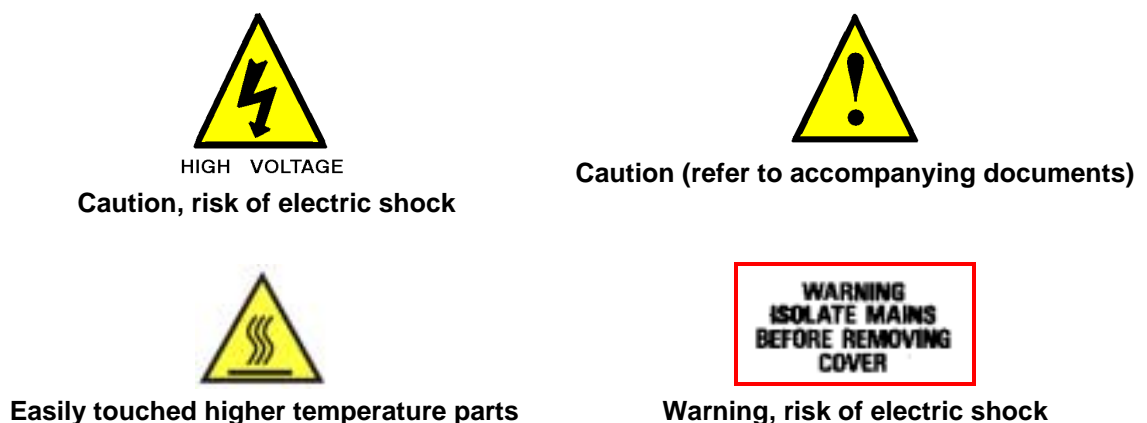


Figure 2.1 - Hazard Warning Symbols

2.5.3 Hazard Warning Labels used in Equipment Manuals

The international warning signs used in equipment manuals as shown in Figure 2.2.



Figure 2.2 - International Warning Symbols

Where appropriate these are used when a specific identifiable risk is involved in either using or maintaining the instrument. These take the form of warning triangles or signs with a graphical description of the hazard.

2.5.4 Instrument Functionality Signs

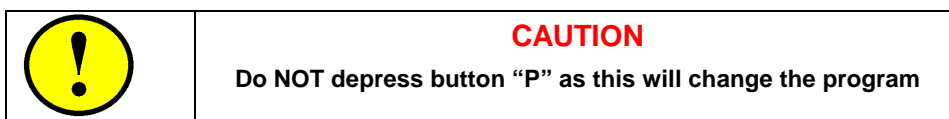


Figure 2.3 - Typical Warning sign as shown in this Manual

This typical sign applies to cautions where there is a risk to the functionality of equipment due to incorrect operation. These cautions or warnings will be contained in a box and be accompanied by a circular warning symbol as shown in Figure 2.3.

2.5.5 Serious Damage to Instruments

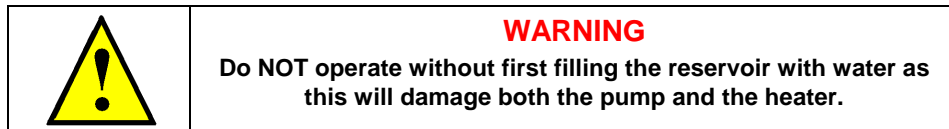


Figure 2.4 - Typical Warning sign as shown in this Manual

This typical caution sign is used where serious damage will be caused by incorrect operation of instrumentation. They will follow the same form as functionality warnings but with a triangular warning symbol as shown in Figure 2.4.

2.5.6 Hazard to Operator



Figure 2.5 - Typical Warning as shown in this Manual

These warnings will generally occur in relevant installation and maintaining sections where there exists a potential hazard to the engineer working on the instrument. They will take the form of the triangular warning symbol accompanied by an international warning sign and bold type lettering beginning with **“WARNING-HAZARD TO HEALTH!”** as shown in Figure 2.5.

2.6 Risk Analysis

2.6.1 Personal Operational Risks

The following is a list of tasks carried out by both the operator and service engineer where recognised risks have been observed, listed is the personnel protection equipment (PPE) which is suggested for use for various tasks on any surface analysis equipment and systems:

Task	Carried out by	Nature of Hazard	Recommended PPE
Cleaning of parts / samples with isopropanol (IPA)	Operator / Service engineer	Splash hazard to eyes, drying of skin	Protective goggles, protective gloves.
Use of Liquid Nitrogen in sample cooling etc.	Operator / Service engineer	Burn risk	Thermally protective gloves and goggles should be worn.
Lifting of Heavy Items	Service engineer	Dropping on foot.	Protective footwear.

Table 2.1 - Personal Operational Risks

2.6.2 Hazardous Materials

◆ Isopropanol (IPA)

For certain service tasks isopropanol is suggested for cleaning components before use in the vacuum system. It should be noted that isopropanol is a flammable liquid and as such should not be used on hot surfaces. In addition it is recommended that protective gloves are worn when using isopropanol.

◆ Compressed Air

Compressed air can be a potential hazard if handled inappropriately. A compressed air line may be fed from some instruments to the customers supply, and the customer should ensure that this and any other service pipes and cables are maintained in good condition.

◆ Nitrogen, Argon and Helium Gas Supplies

Instruments may use nitrogen, argon or helium gas supplies for their operation, the customer is responsible for maintaining the supply to the instrument. This supply should be regulated and kept to the lowest pressure and flow rate as is practical to minimise the effects of any leaks.

◆ Hazardous Gases

Quorum Technologies has no control over the gases used within the system. It is therefore viewed as the customers responsibility to assess the hazards involved and take appropriate precautions when using explosive, toxic or corrosive gases or gases which may result in hazardous products as a result of a chemical reaction.

2.7 Good Working Practices

It is essential that good hygienic working practices are adopted at all times especially in an ultra high vacuum or cleanroom environment and are generally of the "Common sense" type. Some simple good practice rules are:

- ◆ If in doubt don't.
- ◆ If in doubt ask.
- ◆ When handling solvents wear face mask, gloves, apron and work only in a well ventilated area.
- ◆ Mop up any spillages immediately.
- ◆ When handling or decanting mineral oils wear protective clothing.
- ◆ Aerosols of mineral oils, such as that produced by gas ballasting, can prove to be hazardous and an exhaust is recommended.
- ◆ Before attempting to service electrical apparatus, isolate from the mains.
- ◆ Treat all unknown substances as hazardous.
- ◆ Dispose of substances in an appropriate manner.
- ◆ Use the correct tool for the job.
- ◆ Keep a straight back and bend from the knees when lifting heavy objects.
- ◆ Wear protective clothing when using liquid nitrogen.
- ◆ Affix pressurised gas cylinders firmly to walls or racks. Use the correct regulating valves on gas cylinders and always transport cylinders using the appropriate specialist trolley.
- ◆ Obey safety regulations regarding lifts, hoists and machine tools.
- ◆ Always make sure you understand a procedure well before attempting it for the first time.

2.8 SC7610 Specific Safety Hazards

The following Safety Hazards are specific to the SC7610 Sputter Coater.

	<p>WARNING</p> <p>HAZARD TO HEALTH!</p> <p>The Power Supply used in the Model SC7610 unit can operate at up to 1200V D.C</p> <p>HAZARDOUS VOLTAGE OUTPUTS of up to 1200V. D.C.</p>	 <p>HIGH VOLTAGE</p>
---	---	---

2.8.1 Contamination

Contamination can seriously affect the sputtering process. To reduce the possibility of contamination by airborne particles, minimise the time the vacuum chamber is open to the atmosphere.

3 Introduction

This manual is intended for all users of the **SC7610 Sputter Coater** manufactured by Quorum Technologies from the **POLARON** range and provides information on the installation, operation and maintenance of the instrument.

Please note that the servicing and maintenance procedures should only be carried out by qualified service personnel and it is essential that all users should read the **Health and Safety** section of this manual.

3.1 Return of Goods

If goods are to be returned to Quorum Technologies for repair or servicing the customer should contact their local distributor or the factory direct before shipment. A "Returns Authorisation Number" should be obtained in advance of any shipment. This number is to be clearly marked on the outside of the shipment. Complete the returned equipment report form, number **SP-106** with as much detail as possible and return with the goods.

All returned goods are to be accompanied by a completed "Returned Goods Health and Safety Clearance" form **SP-100** attached to the outside of the package (to be accessible without opening the package) and a copy of the forms should be faxed in advance to the factory.

When goods are to be returned under warranty refer to the "Warranty Claim, Repair and Returns Procedure" form number **SP-105**

Copies of all these three forms can be found in the documentation pack supplied with the instrument or direct from Quorum Technologies, the details can be found on page two of this document.

3.2 Return Procedure

Warranty Claim

Electronic and basic servicing capabilities exist at most in-country Thermo companies and at appointed agents, however all components are sold with a **return to factory warranty** (unless otherwise stated) which covers failure during the first 12 months after delivery.

Returns must be sent carriage paid, Quorum Technologies will cover the return carriage costs. This covers defects which arise as a result of a failure in design or manufacturing. It is a condition of warranty that equipment must be used in accordance with the manufacturers instructions and not have been subjected to misuse. This warranty does not cover consumable items such as sputter coating targets and carbon evaporation material. To make a claim under the terms of this warranty provision contact the Customer Service Department at your local Quorum Technologies Representative in the first instance.

Chargeable Repairs

Always contact your in-country Quorum Technologies Representative in the first instance. They will be pleased to assist you and will be able to provide an estimate of repair costs, many offer local repair facilities.

For routine repairs where down-time is not critical. The target standard return time at Quorum Technologies is 20 working days.

Returns

All returns to Quorum Technologies require the following procedure to be followed:

1. Contact the local Quorum Technologies Representative and request a Returns Authorisation Number.
2. Complete a Returned Goods Health and Safety form and returned equipment fault report form.
3. Attach a copy of the completed form to the outside of the package with the usual shipping documents.

Packaging and Carriage

All goods shipped to the factory **must** be sealed inside a clean plastic bag and packed in a suitable carton. If the original packaging is not available Quorum Technologies should be contacted for advice. Quorum Technologies will not be responsible for damage resulting from inadequate returns packaging or contamination of delicate structures by stray particles under any circumstances. All non-warranty goods returned to the factory must be sent carriage pre-paid, (Free Domicile). They will be returned carriage forward (Ex-Works).

4 Description

4.1 Equipment

Each SC7610 Sputter Coater when supplied as a complete package, includes the basic unit, a Gold Palladium target, a start up kit and operation manual. Items can be ordered as a full package or separately against the following numbers:

SC7610 Sputter Coater 220-240V complete, consisting of the following:

LA761002D	Sputter Coater 220-240V. With Gold/Palladium target
SC7610-STARTUP	Start up Kit
OM-SC7610	Operation Manual

SC7610/110V Sputter Coater 110-120V complete, consisting of the following:

LA761003D	Sputter Coater 110-120V. With Gold/Palladium target
SC7610-STARTUP	Start up Kit
OM-SC7610	Operation Manual

4.1.1 Optional items

The following optional items are available from Quorum Technologies:

SC502-314A	Gold Target
SC502-314B	Gold/Palladium Target (1 off supplied with Sputter Coater)
SC502-314C	Platinum Target
SC502-314D	Nickel Target
SC502-314E	Silver Target
SC502-314G	Palladium Target

4.1.2 Accessories

The following accessories are available from Quorum Technologies:

E5005G	Rotary Pump, 90 l/m, 110/240 Volt, 50/60Hz, fitted with E5004 rotary pump exhaust filter.
WS7608	A kit of parts to add a water cooled specimen stage to a SC7610 Sputter Coater, the kit contains: Water Cooled Stage, Inlet and Outlet Connectors, Piping and Fixings.
CA7615	A carbon coating accessory which can be associated with any of the following sputter coaters manufactured by Quorum Technologies is available: SC7620 or SC7640. This accessory which makes use of the vacuum pump and vacuum control facilities of the associated sputter coaters, utilises either carbon fibre or carbon rods in the evaporation process. A 110 volt version is available as part number CA7615/110V
CA076R	Carbon Rod Head for use with the CA7615.
CA076F	Carbon Fibre Head for use with the CA7615.

For further information about the CA7615 carbon coating accessory refer to operating manual OM-CA7615.

4.2 Overview

The **SC7610 Sputter Coater** is a compact magnetron sputter coater, primarily designed for depositing thin conductive metal coatings on to Scanning Electron Microscopy (SEM) samples. Thickness of metal deposition will be determined by the operator, but will typically be in the region of 1 - 20 nanometers.

The SC7610 uses a basic magnetron sputter head with a simple to replace disc target (gold/palladium is supplied as standard, but others metals are available as options). The head is hinged for easy operation and fitted with electrical safety interlocks. The plasma current is variable by adjustment of the vacuum level using an Argon leak valve; the plasma voltage is pre-set.

The 150mm (6") diameter Pyrex cylinder is mounted on an aluminium collar and sealed "L" shaped seals. The small vacuum chamber means pump down times and cycle times are fast; it also allows a small economical rotary to be used. The sample stage is height adjustable over a large range and can easily be removed to accommodate large samples.

For SEM X-ray microanalysis applications the SC7610 can be simply converted to deposit carbon by the addition of an optional carbon evaporation attachment, consisting of a switchable Voltage power supply (CA7615) and a Carbon fibre head (CA076F) or Carbon rod head (CA076R).

The SC7610 comes complete with a one metre of 20mm bore vacuum hose and fittings and requires only the addition of a rotary pump with a capacity of 50 litres / minute or greater (see "options and accessories").

4.3 Technical Specification

4.3.1 SC7610 Sputter Coater Specification

Unit dimensions:	340mm wide x 330mm deep x 330mm high (including vacuum chamber).
Vacuum chamber:	150mm internal diameter x 135mm high.
Weight:	25 kg (55lbs).
Power requirement:	Available for either 230V (13 amp) <i>or</i> 110V (20 amp) operation at 50/60Hz.
Target distance:	Normally 45mm (adjustable).
Power supply output:	Normal operation is up to 800V D.C. at 20mA. Maximum output 1200V D.C.
System control:	Manually by a 180 second timer with 15 second resolution.
Pumping requirements:	Pump to evacuate $>10^{-3}$ mbar.
Sputtering rates:	Refer to Table 6.1, example with a gold/palladium target and current of 25mA a rate of 6nm per minute is achieved.
Coating thickness:	Dependant on time and current, normally between 50 and 300 Angstrom (Å) units for SEM investigations, but will typically be in the region of 1 - 20 nanometers.
Coating uniformity:	Better than 10%.
Gas medium:	Argon.

4.4 Physical Description

The SC7610 Sputter Coater is a simple to operate compact magnetron sputtering system designed to be used for coating Scanning Electron Microscopy (SEM) specimens. The system can only be operated in manual mode.

The SC7610 Sputter Coater is comprised of three main parts:

- The Cabinet assembly
- The Vacuum chamber
- The Sputter head

The cabinet assembly, which supports the vacuum chamber, contains a high voltage power supply, vacuum manifold and gauging. Mounted on the front panel of the cabinet are switches and meters which provide the operator interface to the system. All service connections to the system are made via the rear panel. The cabinet is constructed in a manner so as to comply with the European EMC regulations.

The vacuum chamber is formed by the sputter head top plate assembly, the glass work chamber and the baseplate assembly mounted on the top panel of the cabinet. A safety interlock, conforming to IEC 204, is also incorporated. The integrity of the vacuum is maintained by circular "L" gaskets either end of the vacuum chamber.

The top plate supports the cathode (target) and magnetic deflection system incorporating a dark space shield and a radially magnetised 6 pole magnet which confines the plasma within the target area. Connectors (secured to the top plate), provide the necessary electrical and gas connections between the top plate and the rear panel connectors of the cabinet.

The SC7610 Sputter Coater produces uniform fine grain coating for SEM applications. The sputter head operates at fixed voltage of 800V DC. Coating time is controlled by a 180 second solid state timer with 15 second resolution. The vacuum level and plasma current are monitored by analogue meters. The standard adjustable height specimen stage can be replaced by an optional water cooled stage.

The front panel, includes the chamber pressure and plasma current analogue meters, the timer select switch, the pump start switch, the process (coat) start switch, the set plasma current momentary action switch, and the argon leak and vent valves.

The rear panel incorporates the power on/off switch, the rotary pump electrical outlet socket, the vacuum hose connection port, the argon inlet nipple connected internally to the leak and vent valves, the HV and ground connection sockets, the hose connection nipple for the pressure safety switch and the various fuses. The water connectors are only incorporated when the water cooled stage is incorporated into the baseplate.

For carbon coating (SEM applications only), the top plate assembly can be replaced with an optional carbon rod or carbon fibre evaporation source operated from a low voltage, high current power supply (10V/100A (rod), 20V/50A(fibre)). The sputter head, for safety reasons, is vacuum interlocked to prevent operation of the integral HV Power Supply when the Carbon Evaporation Top plate is used.

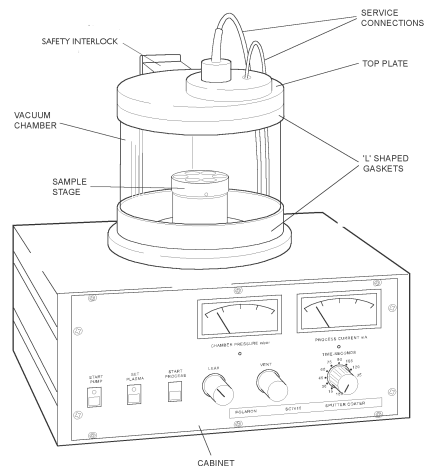


Figure 4.1 SC7610 Sputter Coater

4.4.1 Operational Description

The SC7610 operates at a fixed potential of 800 Volts dc.

An HV voltage is applied between the Target (cathode) and Baseplate (anode) which is at earth potential. A pressure interlock ensures that the HV supply cannot be activated until vacuum chamber pressure is reduced to 10^{-1} mbar or better. Low pressure gas (argon) is leaked into the vacuum chamber to provide a medium for ionisation. Figure 2.2 shows the principles of operation of the SC7610.

Electrons emitted by the cathode, concentrated in the vicinity of the target by the magnetic field, collide with the gas molecules, producing positive ions (due to secondary electron emission). Positive ions attracted by the negative potential of the cathode, bombard the target, causing erosion of the target material. The dislodged target atoms falling toward the sample follow multiple paths due to collisions with the ionized gas, coating the sample on all exposed faces.

A gas discharge glow centered about the cathode is visible.

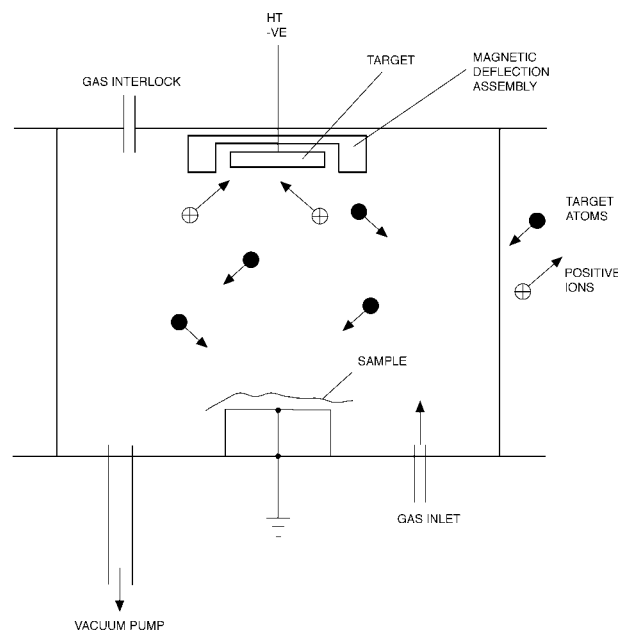


Figure 4.2 Sputter Coater Operation

4.4.2 Pumping Requirements

The work chamber has to be evacuated to 10^{-3} mbar which can be achieved in a reasonable time (depending on the cleanliness of the chamber) using a 90 l/m two stage rotary pump preferably incorporating an anti-suck back device and fitted with an oil mist filter on the exhaust.

4.5 Interlocks

Safety interlocks are incorporated in the SC7610 Sputter Coater to prevent power being switched on with the chamber top plate not in the closed position. Also to cut power to the instrument if vacuum pressure is lost.

	<p>WARNING HAZARD TO HEALTH!</p> <p>Potentially lethal voltages are used in this equipment. Under no circumstances should interlock connections be over-ridden.</p>	 <p>HIGH VOLTAGE</p>
---	---	---

4.6 SC7610 Panel Details

The SC7610 Sputter Coater is designed to be bench mounted, it provides 0 - 1.2 kV D.C. output. Process current of up to 0 - 50 mA is available with a 180 second timer to control the process.

4.6.1 SC7610 Front Panel Controls

The controls and indicators mounted on the SC7610 Sputter Coater front panel are described below and identified in Figure 4.3 and described in Table 4.1.

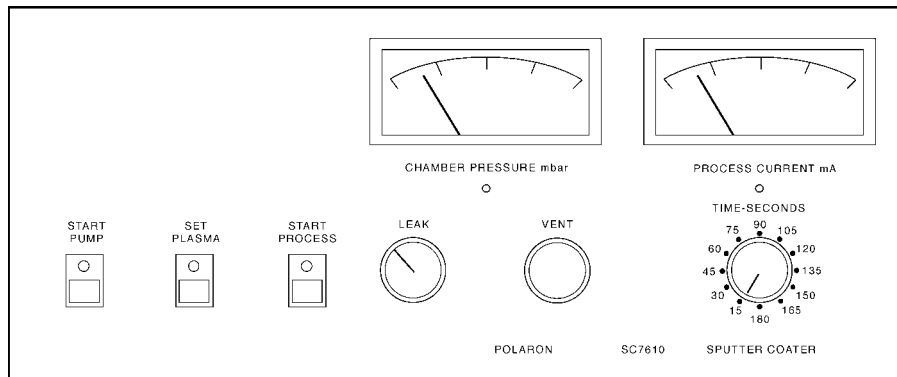


Figure 4.3 – SC7610 Sputter Coater, Front Panel Controls

CONTROL or INDICATOR	DESCRIPTION
START PUMP	A press to operate / press to release switch sets the vacuum pump to ON. After 10-15 seconds a fall in pressure within the chamber will be indicated on the CHAMBER PRESSURE meter.
SET PLASMA	This non-latching switch (hold to operate) is used during the setting of the LEAK valve, which is adjusted to set the correct plasma current.
START PROCESS	Operation of this non-latching switch initiates the sputtering process, the HV supply is activated, until disconnected by the timer.
LEAK	This valve leaks gas into the vacuum chamber to control the plasma current. Used in conjunction with the SET PLASMA switch to adjust the correct operating conditions.
VENT	This valve is opened to admit air into the system after a sputtering process has been completed and the vacuum pump has been turned OFF. During the sputtering process the valve must be SHUT (fully clockwise).
TIME-SECONDS	A twelve position rotary switch sets the required process time, between 15 and 180 seconds, in 15 second steps.
CHAMBER PRESSURE mbar	This meter provides an indication of pressure within the chamber. The normal readings are: Pressure with leak valve closed 0.03 - 0.04 mbar or better. Pressure with leak valve adjusted 0.06 mbar (approx).
PROCESS CURRENT mA	This meter provides an indication of plasma current flow. Under the recommended operating conditions: Plasma current approx. 18 mA.

Table 4.1 – SC7610 Front Panel Control Descriptions

4.6.2 SC7610 Rear Panel Connections

The connection points and other components mounted on the rear panel of the SC7610 Sputter Coater are described below and identified in Figure 4.4 and described in Table 4.2.

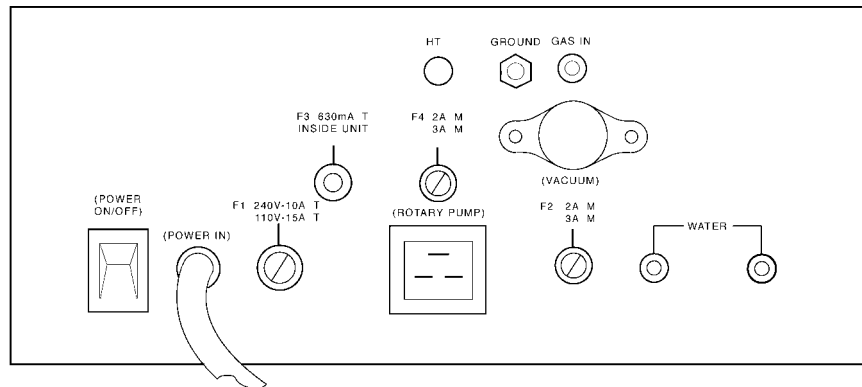


Figure 4.4 – SC7610 Sputter Coater, Rear Panel Connections.

CONNECTION	DESCRIPTION
POWER ON/OFF	A vertically operated paddle switch acts as a mains isolator.
POWER IN	A flying lead provides the connection of power to the system.
F1	Main supply fuse, (240 Volt, 10 Amp, A/S), (110 Volt, 15 Amp, A/S).
F3 (internal)	The 630 mA A/S HT fuse which forms an integral part of the pressure interlock switch. <i>Only accessible when the rear panel is removed.</i>
HT	High voltage socket, connects a negative supply to the cathode mounted in the top plate assembly.
F4	Pirani timer/ controller supply fuse, 2 Amp or 3 Amp.
ROTARY PUMP	A three way mains outlet socket provides power to the rotary pump.
GROUND	Earth connecting point, from the top plate assembly.
GAS IN	A 3 mm nipple provides the gas (argon) input connection point.
VACUUM	A 20 mm vacuum hose connection point, connects to the vacuum pump.
F2	Power supply PCB fuse, 2 Amp or 3 Amp.
WATER	Two 6mm nipples provide Inlet and outlet connections for cooling water supplies. On systems not equipped with water cooling, panel cut-outs are fitted with blind grommets.

Table 4.2 – SC7610 Rear Panel Components

5 Installation

Quorum Technologies has carefully packed the **SC7610 Sputter Coater** instrument so that it will reach its destination in perfect operating order. Do NOT discard any packing materials until the unit has been inspected for any transit damage and the instrument has been used to the customers satisfaction.

If any damage is found, notify the carrier and Quorum Technologies (or local agent) immediately. If it is necessary to return the shipment, use the packaging as supplied and follow the instructions in this manual for return of goods paragraph 3.1.

5.1 Unpacking Checklist

The Equipment package will normally be despatched from the factory in one box. Inside the box the following will be found, refer and check each item off against the supplied packing list.

- ◆ SC7610 Sputter Coater - packed in its own internal packaging, (Target fitted to Top Plate).
- ◆ SC7610 Class Cylinder – packed separately.
- ◆ SC7610 START-UP kit - packed in a polythene bag.
- ◆ Optional Spares - packed individually.
- ◆ Documentation - Inserted in a folder, containing the operating manual and a standard forms pack.

5.1.1 Preparation

- (a) Ensure that a suitable **mains electricity supply** (110 Vac - 20amps or 240 Vac - 13amps, frequency 50/60 Hz) is available. Check that the voltage label attached to the side of the cabinet is suitable for the local voltage and frequency.

The units are supplied for either 230V or 110V operation at 50/60Hz. The power rating is 250VA excluding the rotary pump. The rotary pump outlet is rated at 230V 10A or 110V at 16A. The 240V pump outlet uses a 3-pin plug (404440310) which is supplied or 110V standard US plug (not supplied).

- (b) Ensure that a suitable **gas supply** is available.

Typically: A commercial cylinder of Argon Gas (Zero Grade), fitted with a two stage regulator, in order to deliver gas at a pressure around 5-10 psi (0.7bar).

- (c) Ensure that a suitable **cooling water supply** is available.

This service is not required, unless the water-cooled stage option is fitted.


Ensure that a suitable supply of cooling water is available, a flow rate of 1 - 3 litres / minute of clean water is required. If a closed circuit recirculating chiller is used, avoid operating below dew point.

- (d) Ensure that a suitable **vacuum pump** is available.

Where a rotary pump is used, ensure that the rotary pump has been filled with oil, in accordance with the manufacturers instructions. The exhaust should be filtered or expelled to a safe area. All pumps supplied by Quorum Technologies are fitted with an exhaust filter.

5.2 SC7610 Installation

	<p>WARNING HAZARD TO HEALTH!</p> <p>Potentially lethal voltages are used in this equipment. Before making / breaking connections to the equipment, ensure power is switched off and that it is safe to proceed.</p>	 HIGH VOLTAGE
---	---	---

	<p>WARNING HAZARD TO HEALTH!</p> <p>Precautions to be taken when lifting this equipment. Weight of unit is 25 Kilograms (55lbs)</p>
---	---

- (a) Position the Sputter Coater Cabinet on a suitable level working surface, Access to both front and rear of the cabinet are required.
- (b) Clean the 'L' section gaskets using a cloth moistened with isopropanol and fit to the glass cylinder, then position the cylinder centrally on the base plate (mounted on top of the cabinet).
- (c) Position the top plate assembly on the glass cylinder.
- (d) Position the vacuum pump as close as possible to the Sputter Coater.

5.2.1 Connections

The connections to be made are described below and illustrated in figure 5.1.

- (a) Make the following hose connections to the rear panel of the cabinet.
 - (i) Connect 3mm bore gas tubing (supplied) from argon cylinder regulator to **(GAS IN)** hose nipple.
 - (ii) Connect 20mm bore vacuum tubing (supplied) from the rotary pump to the **(VACUUM)** connector. Ensure the minimum length of hose is used.
 - (iii) If the system is fitted with water cooling make connections to the **WATER** hose nipples using 6 mm bore tubing.
- (b) Connect the Top Plate assembly to the rear panel of the cabinet using the attached connectors:
 - (i) Connect the HV cable, to the **HT** socket on rear panel.
 - (ii) Connect the green/yellow (earth) lead, to the **GROUND** terminal on the rear panel.
 - (iii) Connect the 3 mm bore gas tubing, to the hose nipple (identified as F3) on the rear panel.
- (c) Make the following electrical connections to the rear panel of the cabinet:
 - (i) Connect between the sputter coater power outlet **(ROTARY PUMP)** socket and the vacuum pump.
If the rotary pump is supplied with the Sputter Coater a suitable connecting cable is supplied with the pump, otherwise a suitable 3-way plug (which can be wired by the user) is supplied.
 - (ii) Connect the mains **(POWER IN)** cable via a suitable plug to the local supply, in accordance with the cable colour coding:

Brown	-	Live
Blue	-	Neutral
Green/Yellow	-	Earth


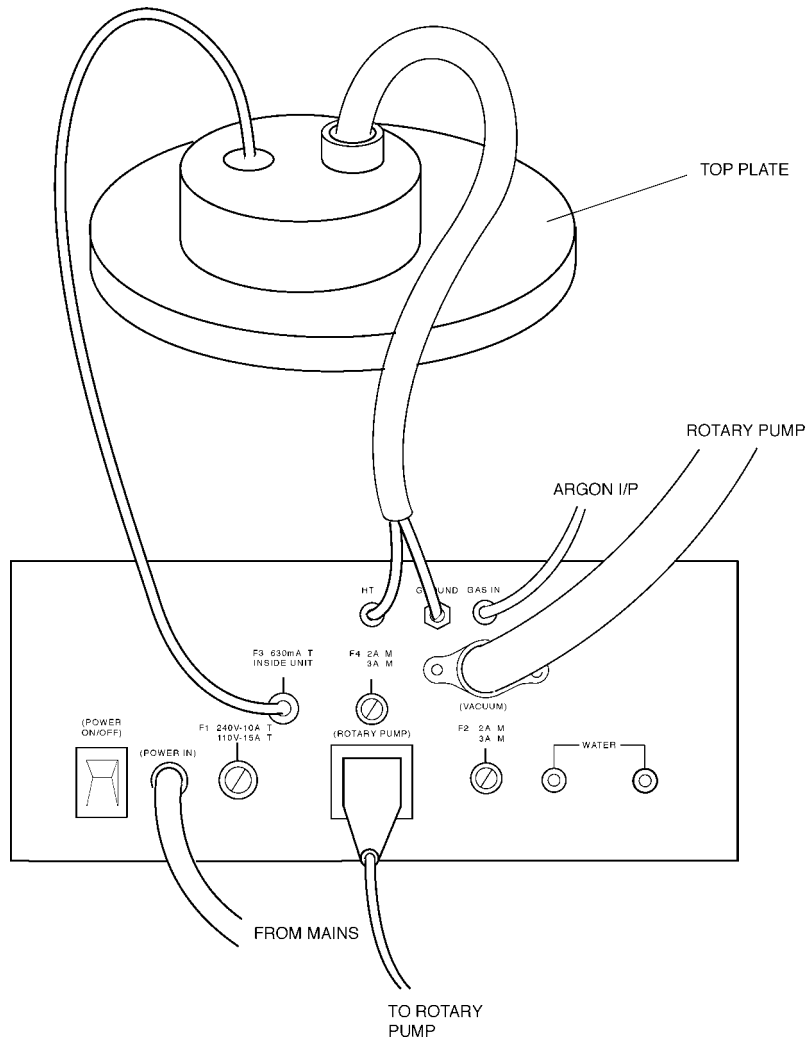


Figure 5.1 – SC7610 Connections

6 Operation

The SC7610 Sputter Coater is designed to be durable and has a long lifetime. The Instrument does contain items, primarily the Target that has finite lifetimes, in order to sustain optimum performance it will need replacing periodically. The lifespan of the Target is dependant on a number of factors, including operating in good vacuum levels, contamination and the purity of the gas being used. The Target should be replaced when it starts to become perforated.

6.1 Test Procedure

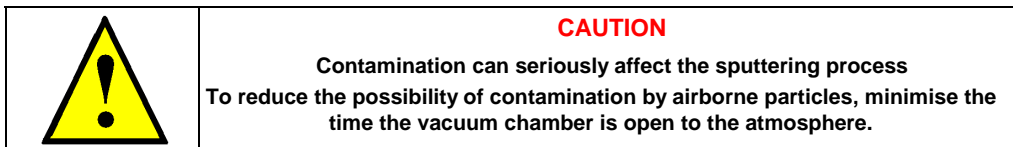

This test procedure, which checks the system is operating correctly, should be performed at the following times:

After the installation process has been completed.

After any operation that could lead to contamination of the vacuum chamber.

(a) Preparation

- (i) Check that **LEAK** and **VENT** valves are closed, (fully clockwise).
- (ii) Check the argon cylinder regulator is open. Set pressure to 0.7 bar (5-10psi).
- (iii) Set **TIME** control to 15 seconds.
- (iv) Check that mains power is available, set the rear panel mounted **POWER ON/OFF** switch to the down position. (ON).

(b) Set vacuum pressure.

- (i) Operate **START PUMP** switch, the integral indicator will illuminate and the rotary pump will start.
- (ii) After 10-15 seconds (dependant on the size of rotary pump), the vacuum gauge will register a fall in pressure within the chamber. Continue pumping until the chamber pressure reaches 3×10^{-2} mbar or better, (indicating there are no major leaks in the system).
- (iii) Open the **LEAK** valve (turn knob counter-clockwise, three to four turns) until the vacuum chamber pressure increases to approximately 6×10^{-2} mbar.

(c) Set process current

Whilst continuing to adjust the **LEAK** valve, intermittently operate the **SET PLASMA** button, the **PROCESS CURRENT** meter will register the current flow and sputtering will occur. Open or close the **LEAK** valve to adjust the current to 18mA. Release the **SET PLASMA** button.

(d) Adjust discharge current

- (i) Operate the **START PROCESS** button, the sputtering process will commence. The current will rise to about 18 mA and a blue or purple discharge glow will be visible, gold-palladium (or other target material) will be sputtered onto the base plate. After 15 seconds, the timer will terminate the process, the discharge will cease and the indicator extinguish.
- (ii) Whilst the sputtering process is in progress, outgassing of the system can alter the process current. Use the **LEAK** valve to correct these changes.

(e) Close down the system, as follows:

- (i) Close the **LEAK** valve (turn fully clockwise).
- (ii) Operate the **START PUMP** button, the indicator will extinguish and the rotary pump will be turned OFF.
- (iii) Open **VENT** valve to admit gas to the vacuum chamber. When sufficient gas has entered the chamber, positive pressure will 'pop' the top plate open
- (iv) Close the **VENT** valve.

6.2 Coating Process

A metal film of uniform thickness between 50 and 300 Å is generally used for SEM investigations.

Care must be taken to ensure the vacuum chamber is kept clean and free from contamination. Contamination which can arise from the out-gassing of specimens, adhesives (especially Chlorohydrocarbon based solvents) and rubber gaskets will adversely affect the quality and rate of sputtering.

A measure of thickness can be obtained using the following equation: **$d = KIVt$**

- d** The coating thickness in Angstrom units.
- K** An experimentally determined constant based on:
 - The metal being sputtered,
 - The gas being used,
 - 45 mm (approx.) target to sample distance.
 - For gold used with argon, K = 0.17 approx.
 - For gold used with air, K = 0.07 approx.
- I** is plasma current, in mA.
- V** is the applied voltage, in kV, (1 kV).
- t** is the sputtering time, in seconds.

For a typical sputtering, using gold in argon with a plasma current of 18 mA for 120 seconds:

$$d = KIVt = 0.17 \times 18 \times 1 \times 120$$

$$= 367 \text{ Å (approx. 3 Å / second)}$$

The uniformity of the coating thickness within the area of the specimen holder is better than 10%.

6.3 Coating Specimens

(a) Mount specimens

- (i) Prepare specimens on stubs, using an approved method.
- (ii) Lift the glass cylinder (with top plate fitted) clear of the cabinet and place aside on a clean surface.
- (iii) Adjust the height of the sample stage. Whilst the most suitable height for a particular application can best be established empirically, 35mm between top of the sample and the target provides a satisfactory general purpose setting.
- (iv) Mount the stubs (with attached samples) on the sample stage.
- (v) Position the glass cylinder and top plate on the sputter coater, to enclose the sample.

(b) Preparation

- (i) Check that **LEAK** and **VENT** valves are closed, (fully clockwise).
- (ii) Check the argon cylinder regulator is open. Set pressure to 0.7 bar (5-10psi).
- (iii) Set **TIME** control to required setting, say 120 seconds (the time required is process dependent).
- (iv) Check that mains power is available, set the rear panel mounted **POWER ON/OFF** switch to the down position (ON).

(c) Set vacuum pressure

- (i) Operate the **START PUMP** button, the rotary pump will start and the reducing pressure will be indicated on the **CHAMBER PRESSURE** meter.
- (ii) When the pressure reaches 0.1 mbar, partially open the **LEAK** valve to flush the vacuum system with Argon. During flushing, allow the pressure to rise to between 0.5 and 1 mbar for 10-15 seconds. Close the **LEAK** valve and allow vacuum chamber pressure to recover to better than 4×10^{-2} mbar.

(d) Set process current

Open the **LEAK** valve and monitor the vacuum pressure. When the pressure begins to rise, intermittently operate the **SET PLASMA** button continuing to adjust the **LEAK** valve. The **PROCESS CURRENT** meter will register the current flow. Open or close the **LEAK** valve to adjust the current to 18mA. Release the **SET PLASMA** button.

(e) Sputter processing

- (i) Operate the **START PROCESS** button, the sputtering process will commence and a blue coloured discharge glow will be visible in chamber. Gold-palladium or other target material will be sputtered onto the samples for the set time period, then the discharge will cease.
- (ii) For a thicker coating, repeat (e (i) above).

(f) Remove Specimens

- (i) Close the **LEAK** valve to shut off the flow of argon to the chamber.
- (ii) Operate the **START PUMP** button, the indicator will extinguish and the rotary pump will be turned **OFF**.
- (iii) Open the **VENT** valve to admit gas to the chamber.
- (iv) When the chamber pressure returns to atmosphere, close the vent valve, lift the top plate clear of the glass chamber.
- (v) Remove the specimens.
- (v) If further specimens are not to be coated immediately, close the top plate to seal the chamber. (See Paragraph 6.3(a)-(v))

6.4 Sputtering Rates

The following sputtering rates have been measured using the FT7690 Film Thickness monitor with a 45mm gap between the targets and the stage crystal holder. The plasma was established using "zero" grade argon gas at the vacuum level required to achieve plasma current. Evacuation of the work chamber was achieved with a 90 l/min two stage rotary pump. Each target was cleaned by sputtering at 25mA for three minutes prior to coating for the measurements.

The graphs show typical sputter rates achieved on a standard instrument under the conditions as specified and are intended only as a guide to coating thickness.


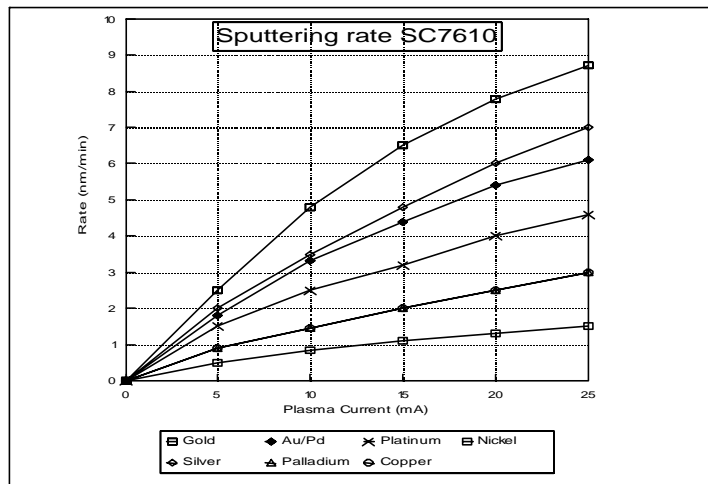


Figure 6.1 – SC7610 Sputtering Rates

7 Maintenance

7.1 Maintenance - General

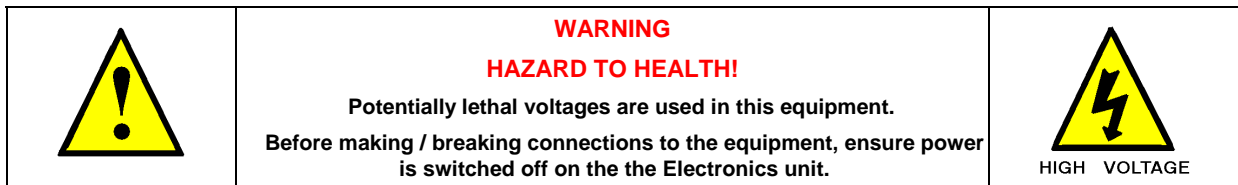

- (a) The procedures listed in this chapter should only be done by persons who have had training and who have achieved a satisfactory knowledge of the necessary skills and techniques.
- (b) If repairs entail the dismantling of any part of the vacuum system, care must be taken to ensure that it is not contaminated (by dust or fingerprints).
 - (i) Always wear disposable plastic gloves.
 - (ii) Do **NOT** handle internal surfaces.
 - (iii) Whenever possible, cover to protect against dust.

7.2 Cleaning

Use a damp cloth or proprietary equipment cleaner to remove surface grime from the outer surfaces of the cabinet and vacuum chamber. Finish with a dry lint free cloth to remove smearing.

7.2.1 Vacuum Chamber Cleaning

Cleaning of the vacuum chamber is required if the interior of the chamber becomes contaminated, cleaning on a regular basis is unnecessary. The fact that the glass walls of the chamber becomes coated with target material and the sample cannot be seen, should not normally be considered reason to clean the system.

If the system is contaminated by handling or air-bourne pollution, carry out the following procedure.

For cleaning use a nylon abrasive pad (Scotchbrite or similar) and Cleaning fluid (Isopropanol or menthyl alcohol).

- (a) Ensure power, water and gas supplies to the sputter coater are set to **OFF**.
- (b) Remove all electrical, gas and vacuum connections from the rear of the instrument. Move the sputter coater to a clean working area.
- (c) Lift the Top Plate assembly clear of the vacuum chamber and place aside. Disassemble the glass cylinder and "L" gaskets from the cabinet assembly.
- (d) Lightly abrade all the accessible interior surfaces of the vacuum chamber to remove any deposits (not the target). Rinse with cleaning fluid. Cover cleaned components to prevent further contamination.
- (e) The target is self cleaning in use and should not require attention. If the target is damaged, replace the target, see section 7.3.
- (f) The vacuum and gas tubing cannot be readily cleaned, if these items are suspect or showing signs of ageing they should be replaced.
- (g) Reassemble the vacuum chamber components.
- (h) Reconnect the electrical and gas tubing connections for the Top Plate assembly.
- (i) Return the sputter coater to its working position. Reconnect the electrical, vacuum, gas and water connectors to the cabinet rear panel.
- (j) When taking the system back into service carry out the Test Procedure (see chapter 6.1), this will ensure the system is thoroughly dried out.

7.3 Target Replacement

In normal use, erosion of the target will occur. Initially this will be seen as a ring on the surface of the target. With further use, holes in the target (generally in a circular pattern) will become apparent. When this occurs the target should be replaced.

7.3.1 Target Removal

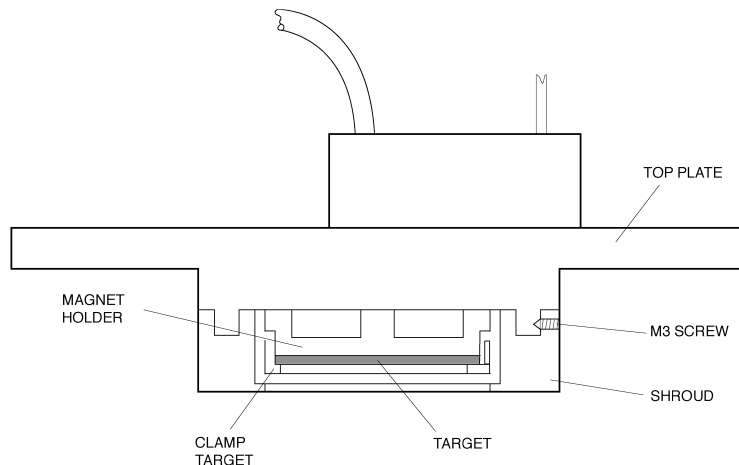


Figure 7.1 – Target Replacement

- (a) Ensure power supplies to the Sputter Coater are set to **OFF**.
- (b) Disconnect the two electrical connections and the vacuum connection to the Top Plate assembly. Cover the open vacuum chamber to guard against contamination by dust. Move the Top Plate assembly to a clean working area.
- (c) Position the Top Plate with target facing upwards, on a clean working surface.
- (d) To release the Shroud, loosen the 3 mm set screw. Remove the Shroud.
- (e) Unscrew and remove the Target Clamp (if the Target Clamp will not unscrew, lubricate lightly with isopropanol). Remove and discard the old target. If necessary clean the Magnet Holder and Target Clamp, using a gentle abrasive material (Scotchbrite), rinse with isopropanol and dry thoroughly.

7.3.2 Refit New Target

For replacement Target type and number, refer to Spares list Paragraph 7.5

- (a) Position the new target on the Magnet Holder, ensure target is lying flat.
- (b) Carefully fit the Clamp over the target, tighten until the target is securely clamped.
- (c) Re-fit the Shroud and secure using the M3 socket screw.
- (d) Position the Top Plate assembly on the vacuum chamber and reconnect the services previously removed.
- (e) When taking the system back into service carry out the Test Procedure (see chapter 6.1), this will ensure the system is thoroughly dried out.

7.4 No Discharge Current

If after switching power on, **SET PLASMA** is operated and the Sputter Coater does not respond, check the following:

- (a) If **SET PLASMA** indicator is illuminated, ensure the vacuum pressure is 10^{-1} mbar or better, if pressure is above this level check for leaks (check **VENT** is closed).
- (a) If **SET PLASMA** indicator is not illuminated check fuse F2 (control supply fuse, fitted to the sputter coater rear panel and F3 (pressure interlock fuse fitted within the sputter coater). If fuse is blown replace with the correct fuse F3, 630mA anti-surge as follows:
 - (i) Ensure power supplies to the Sputter Coater are set to OFF.
 - (ii) Remove all electrical and gas connections to the rear panel.
 - (iii) Remove 10 screws to release the rear panel, retain the screws for future use.
 - (iv) Remove two screws to release the **VACUUM IN** connector, retain the screws for future use.
 - (v) Ease the rear panel away from the main assembly.
 - (vi) F3 fuse holder is mounted on top of the Pressure Interlock Assembly mounted on the rear panel.
 - (vii) Remove the fuse, test and if necessary replace. Discard the blown fuse.
 - (viii) Reassemble the Sputter Coater.

7.5 Spare Parts

Those parts, which due to wear and tear are more commonly required, are listed in the table as follows.

PART NUMBER	DESCRIPTION	QTY
E5004	Oil Mist Filter	1
SC502-314A	Gold Target	1
SC502-314B	Gold/Palladium Target	1
SC502-314C	Platinum Target	1
SC502-314D	Nickel Target	1
SC502-314E	Silver Target	1
SC502-314G	Palladium Target	1
351270380A	Argon Gas Tubing	2 metres
351220320	Vacuum Tubing	1metre
405150310	Fuse – 15A Type T 1 ¹ / ₄ " (F1- 110V)	1
405100310	Fuse – 10A Type T 1 ¹ / ₄ " (F1- 240V)	1
405020210	Fuse – 2A Type T 20mm (F2, F4- 110V)	1
405030210	Fuse – 3A Type T 20mm (F2, F4- 240V)	1
405006210	Fuse – 630mA Type T 20mm (F3)	1

Table 7.1 - Spare Parts

8 Fault Finding

We hope that you experience the minimum of problems throughout the lifespan of the instrument but inevitably problems may occur. Any known problems associated with this type of instrument have been listed below with the possible cause and suggestions to what to do. If problems continue to occur, for example the fuse blows immediately when a new one is fitted, contact the Customer Service Department at Quorum Technologies or your local agent, see Section 9 for list of agents

8.1 Trouble Shooting / Fault Finding

OBSERVATION	POSSIBLE CAUSE	REMEDY	PARA
Unit appears dead	- Cable connections not made	- Check all cable connections	5.2.1
	- Fuse blown	- Replace fuse (F1) on back panel 10A 1 ¹ / ₄ " Anti-surge. - 240V supply <i>or</i> 15A 1 ¹ / ₄ " Anti-surge. - 120V supply	4.6.2
	- Interlock not activating	- Check interlock is operating	-
Rotary Pump not Working	- Unit not switched on	- Check switch on Rotary Pump	-
	- Cable connections not made	- Check all cable connections	5.2.1
	- Fuse blown	- Replace fuse (F4) on back panel 2A 20mm Anti-surge. - 240V supply <i>or</i> 3A 20mm Anti-surge. - 120V supply	4.6.2
Triangular patterns or holes on Target	Target at end of life	- Replace Target	7.3
No Discharge Current	- Low or No Vacuum Pressure	- Leaks in System or Pipework	6.3
		- Vent in Open	4.6.1
	- Fuse blown	- Replace fuse (F3) inside unit 630mA 20mm Anti-surge.	7.4
		- Replace fuse (F2) on back panel 2A 20mm Anti-surge. - 240V supply <i>or</i> 3A 20mm Anti-surge. - 120V supply	4.6.2

Table 8.1 - Trouble Shooting

8.2 Fault Prevention

It is assumed that with a system in regular use and that the system was installed in a suitable environment and in regular use, faults will be repaired as they occur.

To maintain the equipment to the best operating conditions a maintenance schedule is suggested as part of a fault prevention programme, the following items are suggested to be included in such a programme. The frequency of checking will depend on the usage of the equipment.

ITEM	REGULARLY	OCCASIONALLY
Inspect all connections (vacuum, gas and electric) for signs of wear and that they are securely retained in position.	X	
Check all control knobs are secure and operate correctly		X
Check that the lid interlock is operating correctly	X	
Check all earthing cables and connections are secure	X	
Cleaning	X	

Table 8.2 - Fault Prevention

9 Agents

List of Agents Supporting QUORUM TECHNOLOGIES Products.

H D Scientific Supplies Pty Ltd P O Box 6044 Blacktown Business Centre SIDNEY NSW 2148 AUSTRALIA Tel: 61 2962 16955 Fax: 61 2967 12293	Philips Electron Optics BV Building AAE Achtseweg Noord 5 PO Box 218 5600 MD EINDHOVEN HOLLAND Tel: 31 40276 6786 Fax: 31 40276 5141	Medic Corporation Ltd. 20 Peterkin Street Wingate, Private Bag LOWER HUTT NEW ZEALAND Tel: 64 577 0000 Fax: 64 577 2000	Omega Scientific Taiwan Ltd. 5F-1,415 Sec. 4 Hsin Yi Road TAIPEI TAIWAN, ROC Tel: 886 28780 5228 Fax: 886 28780 5225
Labco gmbH Dr - Tritemmel - Gasse 8 A3013 PRESSBAUM AUSTRIA Tel: 43 2233 53838 Fax: 43 2233 53176	OED Technology Limited 22/F Beltrade Commercial Building 3 Burrows Street WANCHAI HONG KONG, CHINA Tel: 852 838 2377 Fax: 852 838 0091	Nerliens Kemisk-Tekniske Aktieselskap Kampengaten 18 N-0654 OSLO NORWAY Tel: 47 2268 5070 Fax: 47 2267 6506	Zen-Yan Scientific Co. Ltd. N° 19, Lane 43 His-Tsun Road Dai-Li City TAICHUNG TAIWAN, ROC Tel: 886 4393 2897 Fax: 886 4393 2401
Soquelec Limited 5757 Cavendish Building Suite 101 MONTREAL QUEBEC H4W 2W8 CANADA Tel: 1 514 482 6427 Fax: 1 514 482 1929	Auro-Science Consulting Ltd Varosfal Kos 5 Postabok 234 1300 BUDAPEST HUNGARY Tel: 361 242 1390 Fax: 361 242 1391	Solid Quest Inc 7 th Floor Westgate Tower Investment Drive Madrigal Business Park Ayala Alabang MUNTINLUPA CITY 1780 PHILIPPINES Tel: 632 850 5759 Fax: 632 850 5961	Becthai Bangkok Equip. Co Ltd. 308/1 Phaholyothin Road Phayathai BANGKOK 10400 THAILAND Tel: 66 2615 2130 Fax: 66 2271 4533
Soquelec Ltd. P O Box 42056 128 Queen Street South Mississauga ONTARIO L5M 4Z0 CANADA Tel: 1 905 569 6613 Fax: 1 905 569 7171	Harley Instruments Plot No. 4, Survey No. 47 Poona Satara Road POONA 411009 INDIA Tel: 91 2042 20602 Fax: 91 2042 20843	Labometer Lda Rue Duque de Palmela No. 30 10 F/G 1250-098 LISBOA PORTUGAL Tel: 351 2135 10180 Fax: 351 2135 25066	Tekser Koll Sti Acibadem Ca. Erdem Sok Baver Apr. No 6/1 Uskudar ISTANBUL 81010 TURKEY Tel: 90 216 327 4041 Fax: 90 216 327 4046
Edrin s.r.o Kon vova 141 130 83 PRAHA 3 CZECH REPUBLIC Tel: 42 2671 08255 Fax: 42 2671 08335	Eisenberg Brothers 13 Gush Etzion St GIVAT SHUMEL 54030 ISRAEL Tel: 972 3532 1715 Fax: 972 3532 5696	Antech Instruments PTE Ltd. 150 Kampong Ampat #05-02 KA Centre SINGAPORE 368324 Tel: 65 289 5500 Fax: 65 289 5522	Agar Scientific 66a Cambridge Road STANSTEAD Essex, CM24 8DA UNITED KINGDOM Tel: 1279 813519 Fax: 1279 815106
Ax Lab A/S Strandboulevard 64 DK-2100 COPENHAGEN DENMARK Tel: 45 3543 1881 Fax: 45 3543 0073	Jeol (Italia) S.p.A. Centro Direzionale Ripamonti Via Dei Tulipani 1 PIEVE EMANUELE (MI) 20090 ITALY Tel: 39 02904 1431 Fax: 39 02904 14303	Set-Point Technology SMM House, 543 Kyalami Boulevard Kyalami Business Park Midrand, 1686 PO Box 11400, VORNA VALLEY SOUTH AFRICA Tel: 27 11466 2200 Fax: 27 11466 2268	ISS Pellowe House Francis Road Withington MANCHESTER, M20 9XP UNITED KINGDOM Tel: 161 445 5446 Fax: 161 445 4914
Elexience ZA Des Godets 9 Rue Des Petits-Ruisseauux 91370 VERRIERS-le-BUISSON FRANCE Tel: 33 1695 38000 Fax: 33 1601 19809	Nanovision Srl Via Treviso, 33 20127 MILANO ITALY Tel: 39 2261 7204 Fax: 39 2261 7215	DAE IL Commerce 2F Jinsung B/D 326-34 Magoek-dong Kangseo-gu SEOUL SOUTH KOREA Tel: 82 2665 8056 Fax: 82 2665 8059	Jeol (UK) Ltd. Silver Court, Watchmead WELWYN GARDEN CITY Herts, AL7 1LT UNITED KINGDOM Tel: 1707 377 117 Fax: 1707 373 254
Gala Instrumente gmbH An der Schmalmach 42 BAD SCHMALMACH D-65307 GERMANY Tel: 49 6124 77952 Fax: 49 6124 60274	Topcon Electron Beam Services Corp 75 - 1 Hasunuma - Cho Itabashi - KU TOKYO 174 JAPAN Tel: 813 3558 2588 Fax: 813 5392 4992	Leica Espana SA Nicaragua 46 BARCELONA 08029 SPAIN Tel: 34 93 49 49530 Fax: 34 93 49 49532	Leo Electron Microscopy Ltd. Clifton Road CAMBRIDGE, CB1 3QH UNITED KINGDOM Tel: 1223 414 166 Fax: 1223 412 776
Asteriadis S.A 56 - 58 Spirou Trikoupi PO Box 26140 GR-100 22 ATHENS GREECE Tel: 30 1 823 5383 Fax: 30 1 823 9567	Hi-Tech Instruments Sdn 60 Jalan Ind. PBP 3 Taman Ind Pusot, Bandar, Puchong 47100 PUCHONG, Selangor MALAYSIA Tel: 60 3589 11638 Fax: 60 3589 11639	Aname C/Jose Fernandex Cuevas 14-B-282244 Pozvelo de Alarcon MADRID SPAIN Tel: 34 91 352 3084 Fax: 34 91 352 3322	Rontec (UK) Ltd Pellowe House Withington MANCHESTER, M20 9XP UNITED KINGDOM Tel: 161 448 0688 Fax: 161 448 0224
Goffin Meyvis P O Box 265, Pergolesilaan 4 4600 AG BERGEN op ZOOM HOLLAND (Also covers, BELGIUM & LUXEMBOURG) Tel: 31 1642 90000 Fax: 31 1642 66651	Prisma Scientific Instruments Services S/B (P.S.I.S.) No. 15B & 15C Jalan Kenari 8 Puchong Jaya Selangor Darul Ehsan PUCHONG 47100 MALAYSIA Tel: 60 3575 2035 Fax: 60 3575 3104	Atema Instrument P O Box 7075 Gillbostraket 8 S-19207 SOLLENTUNA SWEDEN Tel: 46 8626 8365 Fax: 46 8626 8355	Energy Beam Sciences P O Box 468 11 Bowles Road AGAWAM MA01001 UNITED STATES OF AMERICA Tel: 1 413 786 9322 Fax: 1 413 789 2786

10 Index

A ccessories	13	O peration	22
A gents	30	Operators and Service Engineers	8
C hapter Contents	4	Optional Items	13
Cleaning	26	Overview	14
Coating Process	23	P ersonal Operational Risks	10
Coating Specimens	23	Physical Description	15
Conformity	7	Preparation	19
Connections	20	Pumping Requirements	16
Contamination	11	R efit New Target	27
C ontents	3	Return of Goods	12
COSHH	7	Returns Procedure	12
D escription	13	Risk Analysis	10
Disclaimer	8	S afety Policy	7
E quipment	13	SC7610 Front Panel Controls	17
F ault Finding	29	SC7610 Installation	20
Fault Prevention	29	SC7610 Coater Specification	14
G ood Working Practices	11	SC7610 Rear Panel Connections	18
H azard Labels used on Equipment	8	SC7610 Specific Safety Hazards	11
Hazard Signal Words	8	Serious Damage to Instruments	9
Hazard Signals and Signs	8	Service	8
Hazard to Operator	9	Spare Parts	28
Hazardous Materials	10	Sputtering Rates	25
H ealth and Safety	7	T ables	6
I llustrations	6	Target Removal	27
I ndex	31	Target Replacement	27
Interlocks	16	Technical Specification	14
I ntroduction	12	Test Procedure	22
I nstallation	19	Trouble Shooting / Fault Finding	29
Instrument Functionality Signs	9	U npacking Checklist	19
M aintenance - General	26	V acuum Chamber Cleaning	26
M aintenance	26	W arning Labels used in Manuals	9
Manual Layout	3		
N o Discharge Current	28		